

# NON-STERILE GOWNING PROCEDURE

## NON-STERILE GOWNING PROCEDURE BEGINS >



1. GLOVES

Don two pairs of cleanroom gloves, placing the second pair over the first pair.



5. HOOD

Remove hood from packaging making sure that the I.D. label is on the inside. Don hood. Adjust the hood for a comfortable, yet snug fit using the adjusting snaps.



6. COVERALL

Remove coverall from package. Grasp coverall by the neck area to unfold, yet not to touch the ground. Unzip coverall.



9. GLOVES

Roll down donning gloves slightly to expose cuff. Snap cuffs of coverall (if applicable). Place cleanroom working gloves over or under the cuff depending on your requirement.



10. BOOTS

Remove boots from package and don the first boot. Snap upper leg and buckle straps as needed. Swivel that leg to the clean zone without making contact with the top of the bench (if applicable). Don second the second boot. Swivel that leg to the clean zone without making contact with the top of the bench (if applicable).

# NON-STERILE GOWNING PROCEDURE



Don bouffant hair cover and a pair of shoe covers prior to entering the controlled environment gowning area outer corridor.

## 2. HAIR & SHOE COVERS



Enter the change room across tacky mat or other debris control mat taking at least three steps with each foot.

## 3. CHANGE ROOM



Select proper size garments (hood, boots, coverall & goggles)

## 4. SELECT SIZE



Grasp right cuff to bottom right of zipper and left cuff to bottom left of zipper. Gather coveralls upper torso and step into the coverall.

## 7. COVERALL



Tuck in the hood. Completely zip up all zippers. Snap the collar and anklets of coverall (if applicable).

## 8. HOOD



Don goggles. Place goggles so they cover the exposed skin and eye area.

## 11. GOGGLES



Use a mirror to check proper gowning. Discard donning gloves and enter the cleanroom.

## 12. OVERALL CHECK

# NON-STERILE DE-GOWNING PROCEDURE

## NON-STERILE DE-GOWNING PROCEDURE BEGINS >



1. SIT ON BENCH

Sit on the bench. Swivel both legs over to the entry zone without making contact with the top of the bench.



5. REPROCESSING

If you do not intend to re-don the same garment, place the garment in the appropriate receptacle for reprocessing.



6. HANG GARMENTS

If garments are to be re-donned, garments must be hung or stored in a controlled environment.



9. HANG GOGGLES

Hang goggles on hanger or on front snap of coverall. Garment bags may also be used for storage.



10. DISPOSE & EXIT

Exit the change room then dispose of bouffant hair cover, shoe covers appropriately.

# NON-STERILE DE-GOWNING PROCEDURE



Remove goggles. Remove boots one at a time without letting the tops touch the step or floor areas.

## 2. REMOVE GOGGLES



Continue to use a bench to avoid accidental contact with the floor and remove your coverall.

## 3. COVERALL



Grasp cuffs to waist and remove the coverall.

## 4. COVERALL



If coverall is hung snap both boots to the back of your coverall legs so the boot tops do not come in contact with the floor.

## 7. SNAP BOOTS



Remove hood and snap it to the sleeve or neck opening.

## 8. HOOD



End Result. Proper storage configuration for reuse.

## 11. STORE & REUSE

**PLEASE VISIT OUR WEBSITE TO VIEW THE GOWNING AND DE-GOWNING PROCEDURE VIDEO!**

